

# S 3245 UV/VIS DETECTOR

The **Sykam S 3245 UV/Vis Detector** is a variable wavelength UV/Vis detector for routine analysis and sophisticated research. The dual lamp design offers a wavelength range of 190 – 800 nm with a low baseline noise. The front-accessible flowcell can easily be exchanged, as can be the lamps which are accessible through a side panel in the instrument housing.



### ■ Integrated Wavelength Program

The *S 3245 UV/Vis Detector* features a wavelength program to change the selected wavelength over time. With this feature the optimum wavelength can be selected for each analyzed substance according to its retention time.

# ■ Integrated Peak Detector

The integrated Peak Detector works as a basic fraction collector. The peak detection level can be freely programmed for peak start and peak end to enhance the collection purity. An integrated 24V output for switching a solenoid valve is used for the fraction collection, which is automatically operated with a selectable time delay.

## Optional – Dual-Wavelength

The *S 3245 UV/Vis Detector* is available with an optional second wavelength. This feature enhances the Wavelength Program feature that you can measure 2 different wavelengths at the same time. A second D/A converter output comes with this option to keep the system flexible to be used with any data acquisition software available.

### ■ Optional – Online-Scan

Another option for the *S 3245 UV/Vis Detector* is the Online Scan. With the Online Scan whole spectrum information can be gathered at a certain time. This scan information is stored internally and can be accessed at any time. The Online Scan is a good alternative to a full UV PDA detector.



# [mV] 20.0 Peak End Detection Level Peak Start Detection Level Peak Start Time 0.0 1.0 Peak End Time 2.0 Solenoid 3.0 Valve OFF +Delay Fraction Collection Fraction Collection

Figure: S 3245 Peak Detector

Figure: S 3245 Optical Module

# **■ Technical Specifications**

Wetted Materials:	Stainless Steel / PEEK*
Baseline Noise:	± 1 x 10 <sup>-5</sup> AU (@240 nm, 2 sec. Risetime)
Baseline Drift:	< 3 x 10 <sup>-4</sup> AU/h
Wavelength Range:	190 – 800 nm
Wavelength Accuracy:	± 2 nm
Linearity:	> 2.0 AU
Light Source	Deuterium Lamp, Tungsten Lamp
Wavelength Program:	Programmable, 10 steps
Analog Output:	1x 1 V ( <i>optional</i> : 2x 1V)
Control Features:	Internal Peak Detector with +24 V sole- noid switching output.
Dimensions: (W x H x D)	310 x 165 x 478 mm
Power Supply:	100 - 250 ~V (47 - 63 Hz)

<sup>\*</sup> depending on configuration

# Order Information

Catalog No	Description
10 31 011	S 3245 UV/Vis Variable Wavelength Detector
10 31 012	option: Dual Wavelength
10 31 015	option: Online Scan
10 32 001	S 3245 Flowcell, analytical, Stainless Steel
10 32 010	S 3245 Flowcell, analytical, PEEK
10 32 011	S 3245 Flowcell, micro, Stainless Steel
10 32 012	S 3245 Flowcell, micro, PEEK
10 32 013	S 3245 Flowcell, preparative, Stainless Steel
10 32 014	S 3245 Flowcell, preparative, PEEK
40 01 002	Spare Deuterium Lamp
40 01 001	Spare Longlife Tungsten Lamp

### Sykam GmbH

Systeme & Komponenten analytischer Meßtechnik

Gewerbering 15 86922 Eresing Germany

Tel.: ++49 (8193) 93 82 - 0 FAX: ++49 (8193) 93 82 - 20 EMail: Info@sykam.com Web: http://www.sykam.com



